

## QuikRead go Plus Instrument software 2.3.2 release notes

A new QuikRead go Plus Instrument software version 2.3.2 is released on 4.2.2026.

The new software version has been developed and validated in compliance with IEC 62304:2006+AMD1:2015 Medical device software - Software life cycle processes and General Principles of Software Validation; Final Guidance for Industry and FDA Staff, Jan 11, 2002.

The QuikRead go Plus Instrument Instructions for Use (IFU) has been updated with instructions on Bluetooth functionality and other minor improvements. The updated IFU and version history is available on the Aidian [website](#).

### Expanded accessory support

- Added support for Brother QL-800 and Dymo Label Writer 450 Turbo & 550 printers.
- Added support for new barcode readers.
- Added support for Bluetooth-connected accessories:
  - Brother printer models QL-1110NWBc and QL-820NBWC.
  - Bluetooth barcode readers and keyboards.

### Workflow and usability improvements

- Barcode scanners now automatically move focus to the next input field.
- Added a test counter to track the total number of tests performed by the instrument.
- Indicator limits now support decimal values.
- Instrument will reset to admin settings when a general user logs out.
- Updated HbA1c measurement timer to 5.50 minutes.
- Added the option for admin user to delete all QC results.
- Implemented FOB test result USB export in two units.
- Ensured all initial setup parameters are saved.
- Other user interface usability and translation improvements.

**Quality, reliability and connectivity improvements**

- When QC lock is enabled, patient measurements will be locked if QC schedule has expired.
- Prevented using a QC lot that does not match the tested analyte.
- Adjusted rounding in sliding result ranges.
- Improved user notifications.
- Improved error log architecture and overall system reliability.
- Improved LIS and POCT connectivity and WLAN stability.
- Implemented security updates.

**Availability of new software**

After release the new software version 2.3.2 is available on a per customer basis from [www.softwareupdate.quikread.com](http://www.softwareupdate.quikread.com) and is available to order on USB stick (Cat. No 155433).